

# Managing data taken before April 2015

The Portal can also be used to stage data, i.e. copy data from the dCache tape pool to the dCache disc pool. Data can be staged only if the user has a regular DESY account.

During staging, ACLs are created. They grant read-access for the catalogued participants.

- A more detailed view on an archived dataset. This view is obtained by a click on the magnifying glass symbol next to a dataset displayed in the 'Browse data' view. Currently, the dataset may consist of a single or multiple tarballs which can be staged (given a valid DESY account) from here.

**GAMMA-PORTAL**  
Data Management by Physics Groups

Role: USER

Home | Browse data | FTP Registration | Staging status | Migration status | Documentation

Browse data > Files >

Beamtime: 10000307      Period: from 19-NOV-09 to 22-NOV-09

Search   Rows: 15   Actions

select	Archived Directory	Format	Size in Mb	Archived Date ↓	On Disk
<input type="checkbox"/>	space foo	tar	0.74	16-SEP-2013 10:24	no
<input type="checkbox"/>	experiment2_gammadev	tar	0.01	15-FEB-2013 13:18	yes

1 - 2 of 2

Stage   Delete all from disk

\* Stage means to copy multiple datasets to DESY disk space  
\* Delete all from disk means to delete all datasets for beamtime 10000307 from disk (not from tape!)

**Help**  
For downloading, your data must reside on disk. If it is not on disk, please stage it to disk before accessing it. If you do not see the 'stage' button on this page, you are not allowed to use the staging functionality. In this case, please contact the beamline personnel or send e-mail to fs\_ec@desy.de. Use Menu [FTP Registration](#) to select the beamtime you wish to download.

**detailed view**  
-displays users having access to dataset  
-allows staging

- This view is obtained by click on the editing Symbol next to dataset, column 'Users ACLs' from 'Browse Data' view. Depending on the Users role, one see who else has Access to the dataset, who is in Charge of the dataset or one can add/remove users from the list .

**GAMMA-PORTAL**  
Data Management by Physics Groups

Role: USER

Home | Browse data | FTP Registration | Staging status | Migration status | Documentation

Browse data > Users(ACLs) >

Beamtime: 10000472      Period: from 12-OCT-11 to 14-OCT-11

Search: All Text Columns   Go   Actions   Save

	User Role	Door Account	Registry Acco...	Person	Create Date	Change Date	Delete User
<input type="checkbox"/>	applicant	hmustermann	mustermannh	Hans Mustermann	09-FEB-2016 ...	31-JAN-2019 1...	
<input checked="" type="checkbox"/>	leader	testperson	testperson	Test Person	15-NOV-2018 ...	25-FEB-2019 ...	

1 rows selected      Total 2

**Add user to experiment for data access**

Search user   Go

Door account:      Desy account      Pipid

Add User   Add Person for access (visible depending on user role)

**Note**  
To change a role please double-click on the box with the role.

**Help**  
Rights on filesystem:  
**leader** and **applicant**: 'read/write/delete' for all data,  
**participant**: 'read only' to raw directory and 'read/write/delete' for all other data,  
**collaborator**: 'read only' to all data.

### Participant and collaborator

- you will not see the 'Add User' field nor the field / button 'Delete User'. You can not even select roles.

### Applicant

- can add user
- can someone promote to the applicant
- can not downgrade this (or edit)
- can not make someone a leader

### Leader

- can change and delete all roles except 'Leader'.

### Admin

- can change and delete all roles.